

Protocol for Dishwasher in Room 3.7521  
March 2022

The dishwasher is shared across department; sign up on sheet next to S2 autoclave. If someone has signed up after you, call them when finished.

The dishwasher cycle runs in ~50 min.

Alternative racks are available to either side of the room; feel free to carefully rearrange. A setup that often works well for our dishes: flat rack at bottom left for small objects, small bottle/cylinder rack at top left, med/large bottle/cylinder racks middle and right for taller items.

Ensure tall items don't bump into the spinners on roof of dishwasher.

Select "medium" cycle (the name of selected cycle will be flashing) using "select cycle" button, then press "start" twice.

Note that the large (2L+) plastic beakers are a soft plastic that may warp; please wash these by hand. Relatedly—note dishwasher contents may be quite hot when the run is finished.

Lift the door carefully when cycle is done, as contents may have shifted during flight.